Film Degassing Chamber IDEX PN: 9000-2071

Product Datasheet





Description

The 9000-2071 is a 500 µL degassing chamber in a flat film format with ¼-28 flat bottom liquid ports. The fluoropolymer membrane & PEEK chamber is designed to degas solvents at the highest performance levels available for analytical scale HPLC & IC systems. In a small form factor, with reinforced Teflon™ AF film; this degassing chamber has a metal free flow path and universal solvent compatibility, including HFIP and Hexane. The patented chamber design enables low flow restriction when compared to the lumen based chambers.

Degassing Chamber Technical Data

Degassing Channel Membrane	Teflon™ AF, Carbon
Wetted Materials	Teflon™ AF, PEEK, Carbon
Vacuum Housing Material	PEEK
Vacuum Side Materials	PEEK, PFA
Degassing Channel Volume	Nominal 500µl
Vacuum Volume	~100 µl
Maximum Pressure Rating	27 PSIG
Mounting	Threaded 6-32 UNC-2B
Liquid Connections	Threaded 1/4-28 UNF-2B flat bottom, 3/8" depth
Vacuum Connections	Designed for 1/8" ID elastomeric tubing
Operating Temperature	1°C to 60°C (IPA/H ₂ O 20/80)
Storage/Shipping Temperature	-30°C to 70°C
Maximum Flow Rate	10mL/min


Page 1 of 5 2322783 B 12/12/2019






Dimensional Drawing


Dimensions are in Inches (millimeters)

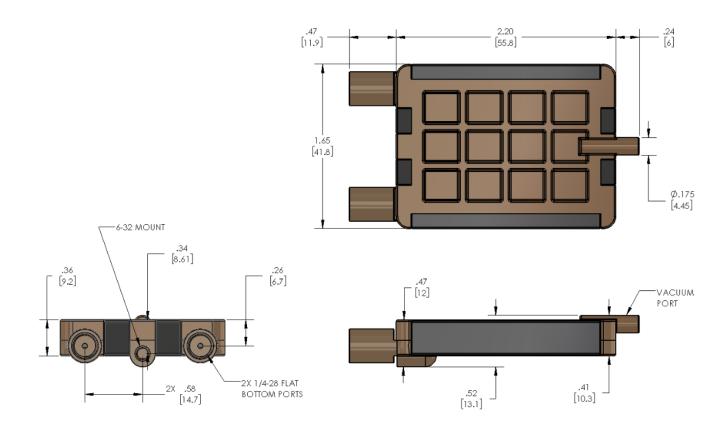

Film Chamber Degassing Efficiency

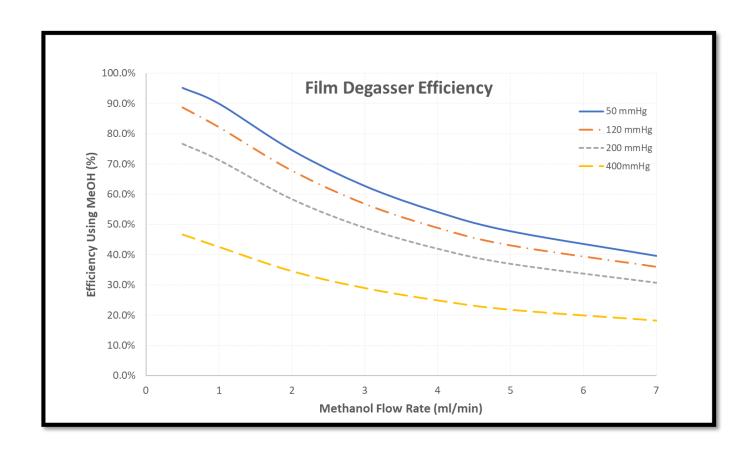

Film Chamber Fluidic Resistance

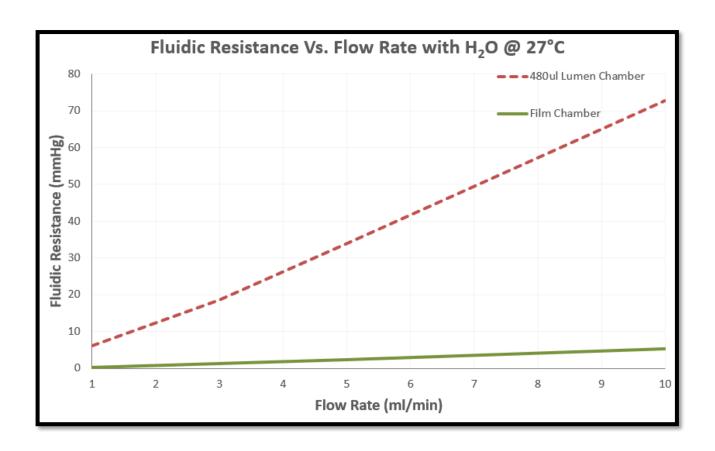

Chemical Compatibility

CHEMICAL COMPATIBILTY		
DESCRIPTION	DEGASSER COMPATIBILITY	
Acetic Acid, Glacial	Compatible	
Acetone	Compatible	
Acetonitrile	Compatible	
Chlorobenzene	Compatible	
Chloroform	Compatible	
Dichloromethane	Compatible	
Dimethyl Sulfoxide	Compatible	
Ether	Compatible	
Ethyl Acetate	Compatible	
Hexafluoroisopropanol	Compatible ³	
Hexanes (60% n-Hexane)	Compatible	
Isopropyl Alcohol	Compatible	
Methyl Alcohol, Anhydrous	Compatible	
Methyl Ethyl Ketone	Compatible	
Methyl Isobutyl Ketone	Compatible	
Methyl tert-Butyl Ether	Compatible	
N,N-Dimethylacetamide	Compatible	
N,N-Dimethylformamide	Compatible	
n-Propyl Alcohol	Compatible	
NMP (N-Methyl-2-pyrrolidinone)	Compatible	
Sodium Azide	Incompatible	
Tetrahydrofuran	Compatible	
Toluene	Compatible	
Water	Compatible	

GENERAL CHEMICALS COMPATIBILTY	
Buffers	Compatible ²
Hydro fluoro solvents	Incompatible
Perfluorinated solvents	Incompatible
Freons	Incompatible
Salts	Compatible ²
Acids	Compatible ¹
Bases	Compatible ¹

Notes:

- 1. Concentrated acids and bases, particularly 1 molar or greater may attack PEEK and PPS wetted parts. Exposure to these acids should be avoided.
- 2. Highly concentrate buffer and salt containing mobile phases should be washed from the HPLC flow path including the degasser flow path if the instrument is to be idle for an extended period. Particular attention should be paid to not leave such concentrated solutions in an unused flow path for an extended period. Crystallization may occur resulting in blockage of the flow path.
- 3. HFIP compatibility is one channel pure HFIP and any number of channels are compatible with HFIP in water.

The above chemical compatibility information has been obtained from our material suppliers, third party sources, and internal evaluations.

IDEX Health & Science LLC has not performed chemical compatibility testing on all solvents listed.

Regulatory Compliance Information

This product is compliant with current RoHS & REACH regulations

Country of Origin

Manufactured & Assembled in USA

Reference Document

Constant Performance HPLC Degassing White Paper Film Degasser Brochure

12/12/2019 2322783 B



