Systec[®] MINI Vacuum Degassing Chamber

The Systec MINI Vacuum Degassing Chamber was designed to be easy to prime, and is configured with a Systec AF[™] degassing membrane to provide maximum degassing capacity with the absolute minimum internal volume (<3% of PTFE designs with comparable degassing capacity).

- Ultra-high degassing efficiency
- Low volume
- Easy primingt

Single Lumen Design


The Single lumen design ensures consistent degassing by avoiding variable flow issues that can be problematic in multi-lumen designs.

Customization Options

Custom degassing capacities, material choices, and metal free designs are available for OEM applications.


Chemical Compatibility

Excellent chemical compatibility with all common solvents used in HPLC. For GPC/Normal Phase LC use degassing chamber 9000-1184.

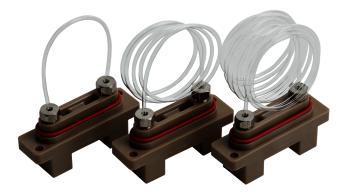



Degassing Efficiency: Residual air for Systec Mini-chambers (percent dissolved air vs. flow rate).

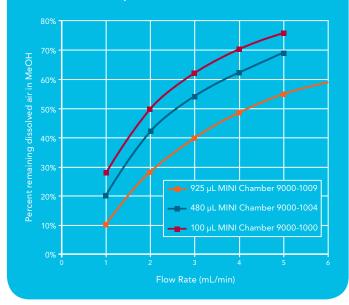

Plot shows remaining dissolved air in methanol using a selection of Systec Mini-Chambers*. The range of chambers and specifications offered provide ample solutions for system designs.

* Water and Methanol mixtures between 30% and 70% methanol will outgas when more than 38% dissolved air remains in each of the solvents. Other water and organic mobile phases being mixed using a low pressure gradient system will undergo similar outgassing.







General Specifications Overall Dimensions Degassing Channel Tubing: Systec AF¹ .50 REF 2.5 REF **Degassing Channel Pressure Rating:** [63mm] [13mm] ▶ 70 PSIG (testing pressure) 0 Wetted Materials: ▶ Systec AF, PEEK, and Glass-filled PTFE or Glass-filled PPS .94 1.75 REF and FEP depending on application [24mm] [44mm] Vacuum Housing Material: Glass-filled PPS (Polyphenylene Sulfide) Hardware Material: ► Stainless Steel or PPS 2.88 **Liquid Connection:** [73mm] ▶ 1/4-28 UNF threaded flat-bottom port Vacuum Connection: Barb-type fitting for 1.57mm (0.062") ID tubing • 自自 00 **ZHCR®** Vacuum Vacuum Pump Controller DEGASSING Dissolved gases are actively CHAMBER removed from a flowing Systec AF™ liquid stream by vacuum via Injector / the Systec AF membrane. Auto Sampler Dissolved gas Fluidic Eluent Pump

Available Configurations^{A,B}

Systec Part Number	Application	Channel Volume (μL)	Max HPLC Gradient Flow Capability ^c (mL/min)	Pressure Drop ^D (kPa/mL/min)	Degassing Flow Path ID (mm/in)
9000-1000	Capillary	100	0.4	0.04	1.14/0.045
9000-1004	Analytical	480	2.0	0.18	1.14/0.045
9000-1006	Analytical	670	2.8	0.25	1.14/0.045
9000-1009	Analytical	925	4.0	0.36	1.14/0.045
9000-1184	Analytical/GPC	480	2.0	0.18	1.14/0.045
9000-1810	Analytical/GPC	285	1.5	No data	0.89/0.035
9000-1818	GPC	925	4.0	0.36	1.14/0.045

A. Custom configurations are available. Consult the factory for your specific application.

B. The standard MINI degassing chamber is not recommended for GPC applications (room temperature or heated), nor for use with HFIP (hexafluoroisopropanol). We offer a special GPC "hardened" version. Please consult the factory for details.

C. The flow rates given are for a gradient mixture of 50/50 MeOH/H2O, with a typical low pressure gradient mixing valve. Higher flow rates are possible with high pressure mixing. D. Estimated tubing pressure per unit change in flow assuming laminar flow with a viscosity of 1.0 cP.

HEALTH & SCIENCE



Global Premium Distributor: Biotech AB - Europe | Biotech LLC - USA | Bionik Inc - Japan www.biotechfluidics.com info@biotechfluidics.com | +46 300 56 91 80